




Phenmetrazine

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.

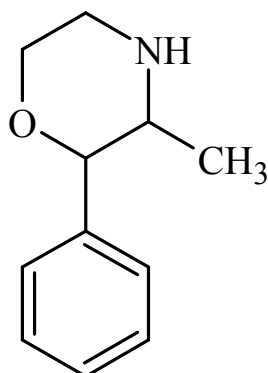

1. GENERAL INFORMATION

IUPAC Name:	3-methyl-2-phenylmorpholine
CAS#:	134-49-6
Synonyms:	Not Available
Source:	DEA Reference Material Collection
Appearance:	White powder
UV_{max}(nm):	Not Determined

2. CHEMICAL AND PHYSICAL DATA

2.1 CHEMICAL DATA

Form	Chemical Formula	Molecular Weight	Melting Point (°C)
Base	C ₁₁ H ₁₅ NO	177.24	Not Determined
HCl	C ₁₁ H ₁₅ NO HCl	213.70	173.8



Phenmetrazine

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.




3. QUALITATIVE DATA

3.1 NUCLEAR MAGNETIC RESONANCE

Sample Preparation: Dilute analyte to ~23 mg/mL in D₂O containing TSP for 0 ppm reference and maleic acid as quantitative internal standard.

Instrument: 400 MHz NMR spectrometer
Parameters: Spectral width: at least containing -3 ppm through 13 ppm
Pulse angle: 90°
Delay between pulses: 45 seconds
1HNMR: Phenmetrazine HCl; Lot# 08/10/13, D₂O; 400MHz

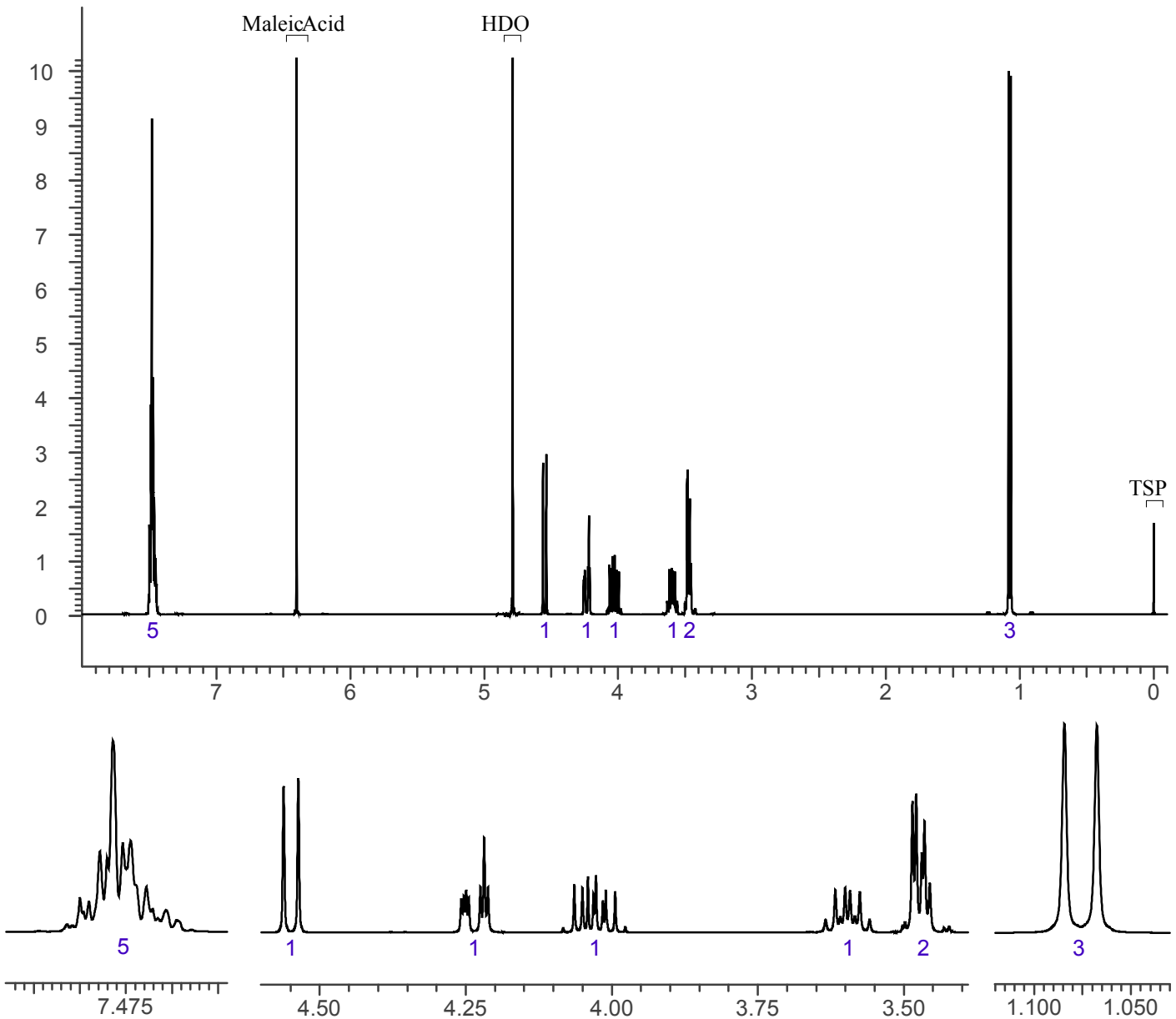



Phenmetrazine

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.



3.2 GAS CHROMATOGRAPHY/MASS SPECTROMETRY

Sample Preparation: Dilute analyte ~4 mg/mL in CHCl₃ base extracted with sodium carbonate.

Instrument: Agilent gas chromatograph operated in split mode with MS detector

Column: HP-5 MS (or equivalent); 30m x 0.25 mm x 0.25 μm

Carrier Gas: Helium at 1.5 mL/min

Temperatures: Injector: 280°C

MSD transfer line: 280°C

MS Source: 230°C

MS Quad: 150°C

Oven program:

1) 100°C initial temperature for 1.0 min

2) Ramp to 280°C at 12 °C/min

3) Hold final temperature for 9.0 min

Injection Parameters: Split Ratio = 25:1, 1 μL injected

MS Parameters: Mass scan range: 30-550 amu


Threshold: 100

Tune file: stune.u

Acquisition mode: scan

Retention Time: 6.582 min

EI Mass Spectrum: Phenmetrazine HCl; Lot# 08/10/13

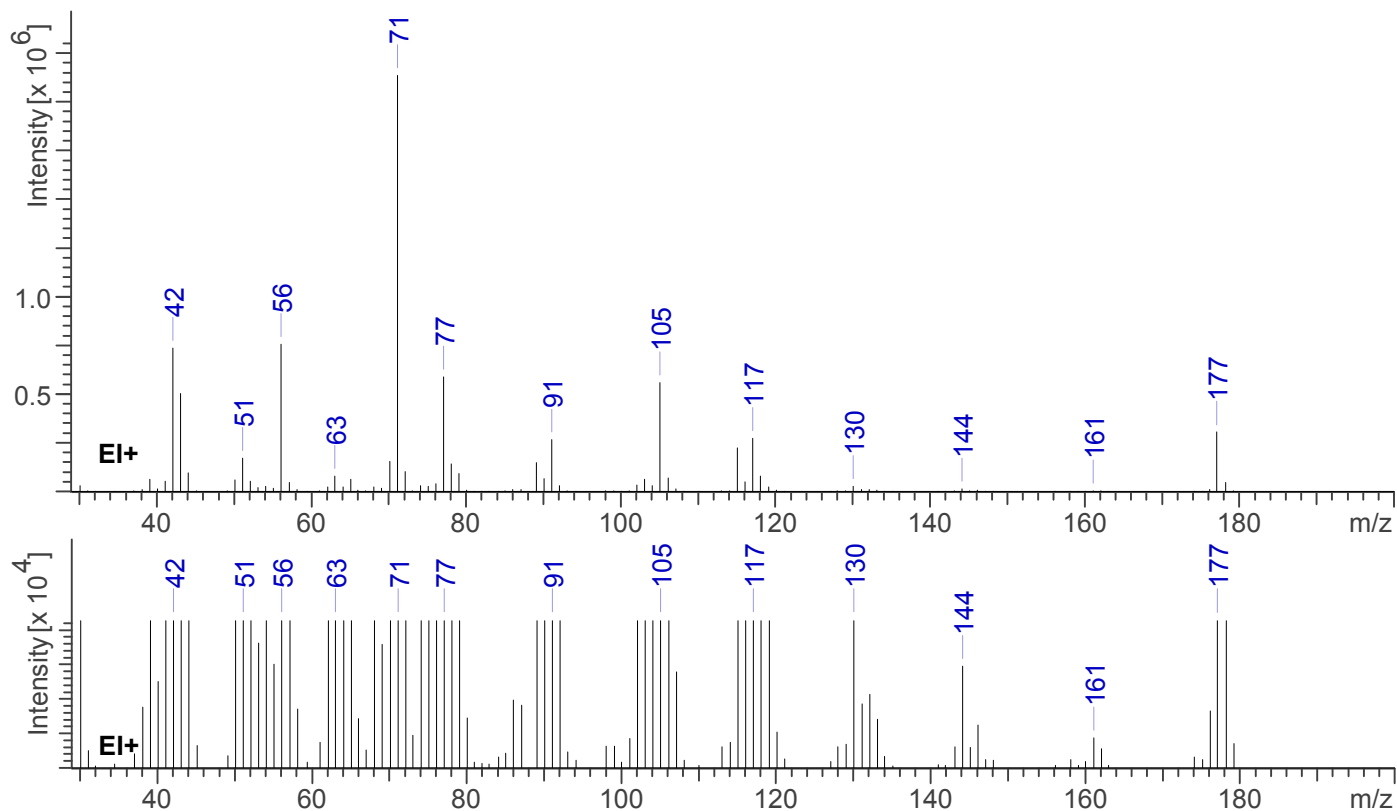



Phenmetrazine

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.





3.3 INFRARED SPECTROSCOPY (FTIR)

Instrument: FTIR with diamond ATR attachment (1 bounce)

Scan Parameters:
Number of scans: 32
Number of background scans: 32
Resolution: 4 cm⁻¹
Sample gain: 8
Aperture: 150

FTIR ATR (Diamond 1 Bounce): Phenmetrazine; Lot# 08/10/13

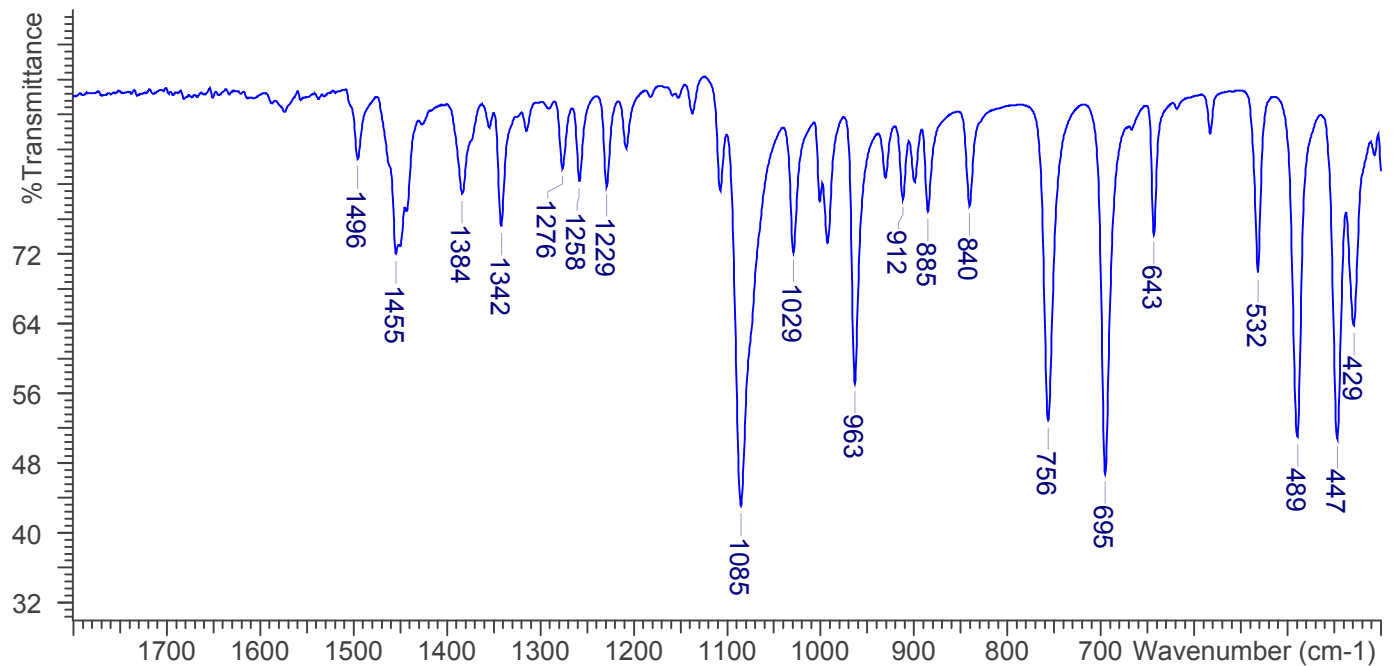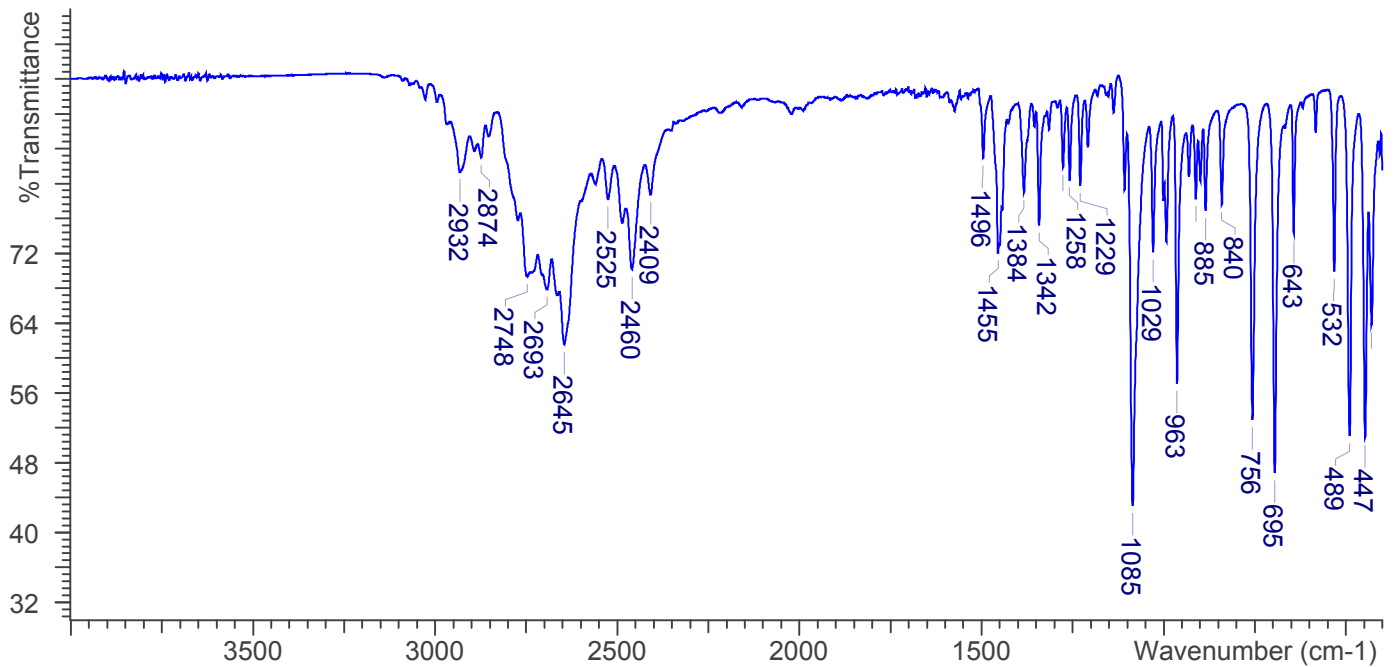



Phenmetrazine

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.



4. ADDITIONAL RESOURCES

[*Forendex*](#)

[*Wikipedia*](#)